# ZOOSCAN

## MANUAL TO SCAN and PROCESS SAMPLES

### Quantitative Imagery Platform of Villefranche sur Mer (PIQv)

Picheral, Elineau

2018/09/25

*J. Plankton Res.* Gorsky et al. 32 (3): 285. : ZOOSCAN methodology paper.

Open access link in [www.zooscan.com](http://www.zooscan.com)

<http://plankt.oxfordjournals.org/cgi/reprint/fbp124?ijkey=gJnvV9AjR20gITg&keytype=ref>

## 1. IMPORTANT NOTE ON ZOOPROCESS VERSION

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## 2. ZOOSCAN WORKING PROTOCOL ZooSCAN 工作手册

This working protocol might be adjusted by users according to their experience and needs in order to take advantage of the tools provided by the application. We anyhow recommend to always keep the default settings offered by Zooprocess.

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The typical analysis sequence is as follow: 典型的分析顺序如下：

- Fill in Sample metadata (as many sample as useful)  
填写样品元数据信息（可用样品越多越好）
- SCAN Background (on a daily base) 扫描背景（在每日基础上）
- SCAN sample (one or several in a raw) 扫描样品（在原始文件夹中一个或几个）
- CONVERT and PROCESS samples 转换和处理样品
- CHECK process 检测流程
- SEPARATION of touching objects 将彼此接触的对象分离开
  - o On vignettes (best option) 在小图上（最好的选择）
  - o On global image, either B&W or in grey level (old option)  
在球型图片上，在 **B&W** 或在灰度值水平上（以前的选择）
- PROCESS image again to include separation mask and get a better final dataset  
再次处理包含 **separation mask** 文件夹在内的图片，以便获得更好的终极数据库
- IMPORT in Ecotaxa to predict identification and manually validate or correct the prediction made by the application.  
导入到 **Ecotaxa** 来进行预测识别，并通过手动验证或更正由应用程序所做的预测。

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- Open Image J. Zooprocess starts (or click on the Z icon if ImageJ is already running) and choose the option “CREATE NEW PROJECT” which is at the bottom of the project list.

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- Choose the drive. 选择一个存储盘
- Select the scanning option: we recommend choosing the Large frame for images at 2400 dpi resolution. The Narrow frame for images at 4800 dpi resolution can be utilized only if you run a ZooScan V3 or V4 with a Windows 7 pro, 8 or 10 64

bits computer. It does not provide much better resolution and limit the scanning area and the number of organisms in the image.

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*A folder will be created on the selected drive and all files from this project will later be saved in the specific subfolders.*

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#### 4. Fill in Sample metadata 填写样品信息

This operation is no longer connected with the scanning of the samples. You can fill in metadata in advance for several samples. You can even do it on the field and then move your project to the computer which is connected to the Zooscan.

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At this point you should have carefully considered the naming convention of your samples! As it not possible to change the name of your sample afterward. And you should carefully check the format of several metadata such as Latitude and Longitude (see the metadata windows)!

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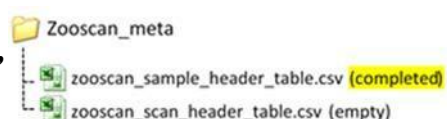
*An excel table named :*

*“zooscan\_sample\_header\_table.csv” will be created by the application in the “Zooscan\_meta” folder of your project. This table lists all the samples and associated metadata you have filled in.*

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You can modify metadata (except the sample name) any time between all the steps clicking the option “EDIT and MODIFY metadata”.

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You can EDIT metadata at any time to correct possible mistakes or complete information on the samples or the scans. Pay attention at these metadata. They will be essential for the later analysis of the results.

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*The tool corrects the metadata in the tables of the meta folder and inside all files related to the sample anywhere in the project.*

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## 6. SCAN (CONVERT) Background Image 扫描（转换）背景图片

The background image is a “blank” image that will be used during the image analysis process. It should be made before the samples are scanned and using the same parameters than for the scan of the samples. You have to scan a minimum of two background images (just follow the steps!) that will be combined in a unique “blank” image. You will do it at the beginning of every scanning session when you turn on your ZooScan, so usually every morning before the first sample.

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*The two background images will be processed into a unique background image and automatically saved in “Zooscan\_back” folder.*

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1. Turn on the ZooScan (biotom, v1 and v2: power button on grey box; v3 or v4: power button behind the instrument), manipulate the ZooScan gently and maintain it (e.g., always hold the tray when you lift it to remove the water or it could fall).  
打开 ZooSCAN 电源（biotom, v1 和 v2 版本：电源按钮在灰色盒子里；v3 和 v4 版本：电源按钮在仪器的后方），轻柔地操作和维护 ZooSCAN（例：当你把样品池举起来倒掉水的时候，一定要一直扶着它，否则它可能会掉下来）。
2. Clean and rinse the scan tray and the cover glass using freshwater to remove the dirtiness.  
用淡水清洗和冲洗扫描池和盖板玻璃，以清除污垢。
3. Eliminate marks on the glass and the frame and check from time to time if the glass of the ZooScan cover has no mark.  
清除玻璃和样品框上的污垢，并不时检查 ZooScan 盖子上的玻璃是否没有污垢。
4. Pour some clean freshwater (stored at room temperature) to cover the tray.  
倒一些干净的淡水（室温保存）盖住样品池。
5. Place the frame that was defined for your project (NARROW/LARGE). Control that the frame is well placed on the recommended and tagged corner of the scanning tray (the area of scan is set up to include the frame borders). For biotom, v1 and v2 instruments there is a mark on the frame for his placement on the tray. Check the manuel delivered with the ZooScan.

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6. Fill the tray with freshwater until the edge of the frame is covered. Check that there is no dust or bubble below the OD (dark circle) position and that the tray and the water are clean : remove dirtiness and bubbles using a plastic pipette or a cactus stick provided with the instrument to avoid scratching the glass tray.

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Figure 1 : Filling the Zooscan tray

7. BIOTOM, V1 and V2 models: DON'T FORGET to turn on the light (green button or lever button) just before scanning. V3 and V4 models: Check that the light rotating switch (left side of the Zooscan) is on UPPER position.

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8. Click on “SCAN (CONVERT) Background Image” tool. Follow the instructions that pop up on your computer screen. Wait the recommended time (30 sec.) between the preview, scan n<sup>o</sup>1 and scan n<sup>o</sup>2 (cf. the instructions window on your computer screen). Don't forget to press OK in the instructions window before the second scan.

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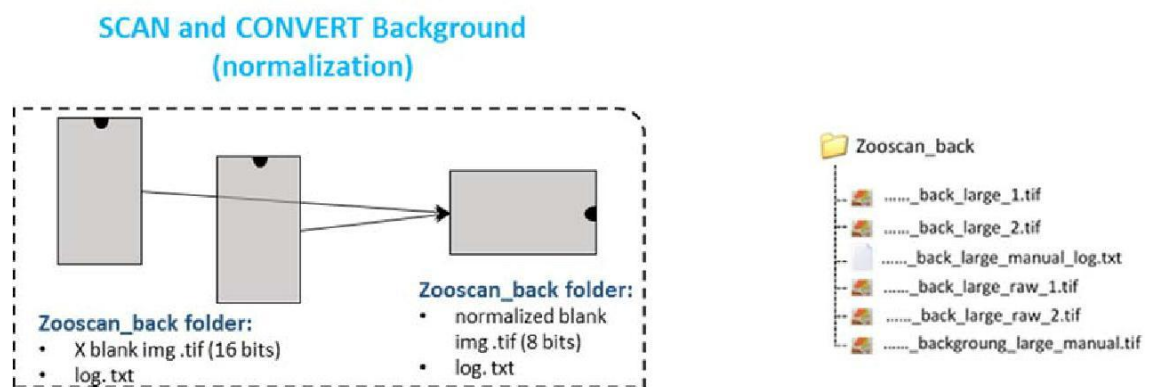
Figure 2 : Process of background images

## 7. Sample preparation and scan 样品准备和扫描

### 7.1 Sample preparation 样品准备

**Please do consider that the FORMALIN is a carcinogen and mutagen product. Apply all necessary safety procedure to remove all risks when you remove the formalin from the sample:**

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**For safety reasons, we never pour directly the samples containing formalin on the Zooscan tray but we previously remove the seawater and formalin from the sample and use filtered seawater (or fresh water) instead.**

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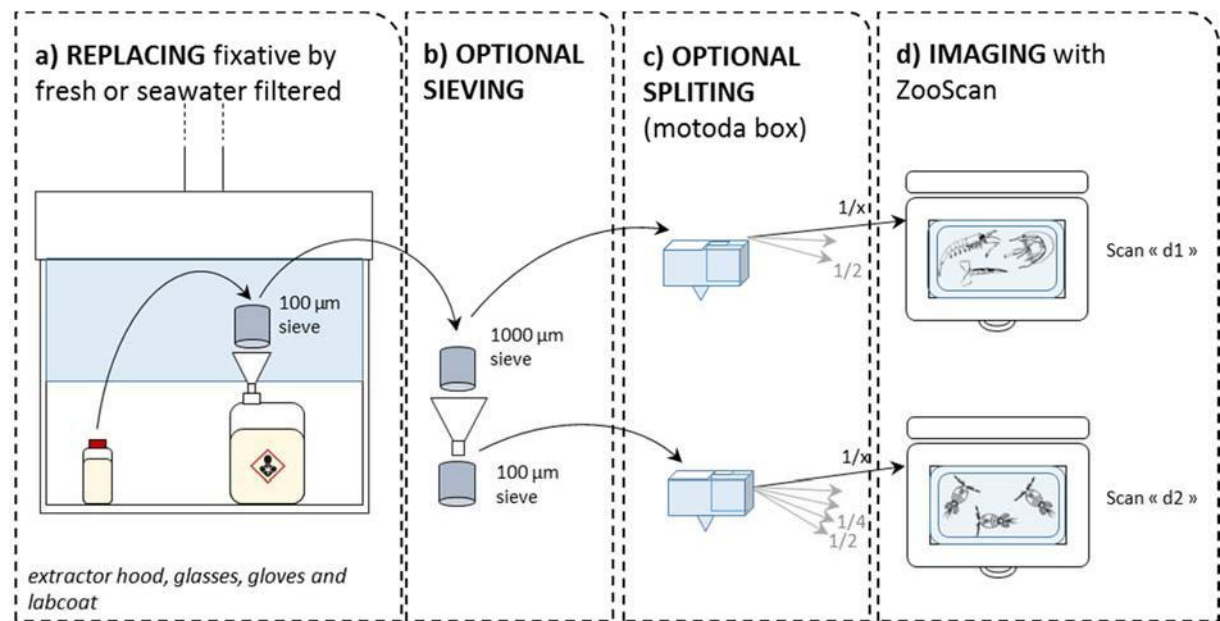
Figure 3: Procedure to prepare a sample collected using a WP2 net (200  $\mu\text{m}$ ): a) removal of the seawater and formalin, b) the sample is optionally sieved in two fractions, c) each of them is optionally split in order to get convenient numbers of organisms and d) scan of both fractions ("d1" and "d2")

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The sample is sieved under the extractor hood to remove the preservative and the sea water.

You can keep it to re-fill the sample afterwards.

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In order to prevent the underestimation of large and rare organisms and depending on the mesh size of the net utilized to collect the sample, the sample is then sieved through a 1000  $\mu\text{m}$  mesh and 100  $\mu\text{m}$  mesh to get 2 size fractions.

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The two fractions are then split before scanning to obtain a proper number of organisms: a large one (>1000  $\mu\text{m}$ ) and a small one (100  $\mu\text{m}$ -1000  $\mu\text{m}$ ).

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**NOTE:** 注意：

- You can control that it is useful to sieve the sample in two size fractions by comparing the splitting ratio necessary to obtain a good image or a correct quantity of organisms for the two fractions. **This ratio must be higher for the lower fraction.**

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- By experience we scan 500-1000 organisms for the large fraction and 1000-1500 organisms for the the lower fraction. These numbers can be adapted according to user experience and care. The number of touching organisms must be as limited as possible to avoid spending much time to later separate them from the images.

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- You can decide to split the whole sample instead of using two size fractions (net mesh size > 500  $\mu\text{m}$  for example).

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- In order to avoid mistakes during the splitting procedure using the Motoda box, we recommend to keep all successive remaining fractions in pre-labelled jars (1/2, 1/4, 1/8, 1/16....).

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## 7.2 STEP BY STEP SCANNING PROCEDURE 一步接一步的扫描过程

### NOTES: 注意:

- The METADATA must be previously filled (see above)  
元数据信息必须提前填好（参见上文）
- The BACKGROUND must be scanned (see above)  
背景图像必须扫描（参见上文）

1. Pour some water (freshwater or filtered seawater) on the scanning tray until the glass is totally covered (don't cover the edges of the frame at this step as the addition of the sample will then cover the edge). In order to prevent condensation and bubbles to appear on the ZooScan tray (because of the different temperatures between the working room and the water in the pipes), we recommend to anticipate your water usage and store few liters of water in the same room of the instrument.

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2. Place the frame. The frame size (NARROW or LARGE) depends on your choices at step 2 "Create new project". Check that the frame is well placed, touching the recommended and tagged corners of the scanning tray as for the scan of the background (see above).

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3. Clean water droplets or marks on the frame and remove the dirtiness in you water with pipette and the cactus sticks.

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4. Pour the sample from the last fraction of the Motoda box  
倒入 **Motoda** 分样器中最后一次稀释的样品。
5. Add water (if necessary) until all the edges of the frame are covered as for the background. If you pour too much water, the floating organisms will be out of focus. You thus have to struggle between : enough water to cover the edges and an insufficient level to avoid out of focus organisms.  
加水（如果必要的话）直到样品框的所有边缘被覆盖（如背景扫描步骤）。如果你倒了太多的水，漂浮的组织将会失去焦点。因此，你必须在两点间取得平衡：有足够的水覆盖边缘，又不至于过多使组织失去焦点。
6. Allow 5-15 minutes to separate touching organisms with the cactus sticks. No organisms on the edge and along the edge! Pay attention to the separation of the objects from each other. Make a compromise between the time spent for this task and the quality of the image. After the process of the image, you can separate touching objects on the final image with the separation tool (see below) in Zooprocess but you may lose details on the organisms and they may be truncated.  
花 5-15 分钟时间用仙人掌刺分离彼此接触的组织。不要有组织在透明框的台阶上或者在台阶附近！在将对象彼此分离的时候要多加小心。花费在这项工作上的时间与图片的质量之间，你要做好取舍。在图片处理之后，你可以用 **Zooprocess** 中的分离工具（见下图）分离最终图片上相互接触的物体，但是你可能会丢失生物的细节，它们可能会被截断。
  - You can accelerate the separation of organisms by pouring the sample homogeneously on the tray and by pouring the extra water on conglomerated organisms of the tray to “dissolve” them.  
你可以通过将样品均匀地倒在样品池上，以及在样品池中聚集生物较多的地方多加一些水来“分解”它们，从而加速样品的分离。
  - Place the larger individuals in the center of the tray because the image will be cropped on the edges of the frame. If some organisms are floating, try to sink them by little pushes with the cactus sticks (the size measurements of floating organisms are biased, and their image captions are blurred). If they do not sink and are very few, the best option is to move them apart of the image (e.g. placing them on the edge of the frame). This step is critical to have good data quality.  
将较大的个体放在样品池的中心位置，因为如果在透明框的边缘，图像将会被裁剪。如果一些动物浮在水面上，试着用仙人掌棒轻压它们，使它们下沉（动物漂浮着的话，尺寸测量会有偏差，并且它们的图片会模糊）。如果它们不下沉的话并且又比较少，那最好的选择就是把它们从图片中分离出来（例如把它们放到样品框的台阶上）。这一步骤对于拥有良好的数据质量来说至关重要。
7. Check that there is no condensation on the glass of the ZooScan cover.  
检查在 **ZooSCAN** 上盖的玻璃上没有冷凝水
8. Launch Zooprocess, select your project and click on the option: “SCAN sample with Zooscan (for archive, no process)”.  
打开 **Zooprocess**，选择你的项目并点击“**SCAN sample with Zooscan (for archive, no process)**”选项。
9. Select the sample name. Then, follow the instructions that pop up on the screen (Zooprocess Instructions window). In Zooprocess, select the proper fraction name “d1”, “d2”, “tot” or user defined according to the fraction ( $100\mu\text{m} < \text{d}2 < 1000\mu\text{m}$ ,  $\text{d}1 > 1000\mu\text{m}$ , “tot” if no sieving) when asked.

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10. Follow carefully the instructions in the Zooprocess windows to start the scan of the sample (one scan only). Do not forget to press OK in Zooprocess after pressing SCAN in Vuescan. Do not touch any key of the keyboard during the scan ! Do not make any vibration on the Zooscan that may perturbate the water surface on the tray! Biotom, v1 and v2 Zooscan versions : do not forget to turn on the green light of the Zooscan before launching the scan in Vuescan. WAIT 30 seconds between the preview and the actual scan. v3: select the UPPER light!

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#### ADDITIONAL INFORMATION: 附加信息:

- Three files will be created in the “\_raw” folder of the project. The filename is composed of the sample name plus the fraction name (sample\_d1.\*) 三个文件将被创建在项目的“\_raw”文件夹下。文件名字由样品名称加上分级名称 (sample\_d1.\*) 组成

- TIF image (16 bits) will be processed later when using “CONVERT & PROCESS Images and organisms in batch mode” option in the Zooprocess main menu.

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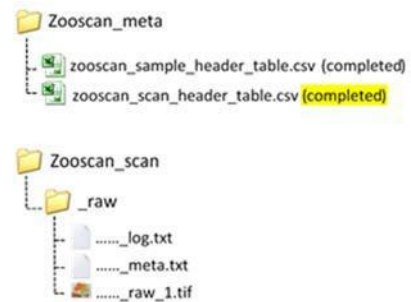
- META file gives information on the sampling method (e.g., sampling site net dimensions, tow, volume) and on the sample preparation for the Zooscan (e.g., pre-filtering and subsampling ratio). It summarizes the information contained in the sample and scan tables from the meta folder of the project.

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- LOG file records information on the scanning method (parameters).

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## 8. RECOVER sample 回收样品

1. Remove and rinse the transparent frame with a squeeze bottle above the scanning tray to recover all specimens that may be stuck on it.  
取出样品框，并使用能挤压的洗瓶在样品池上方清洗样品框，以回收所有可能粘在其上的样品。
2. Remove the sample from the tray: lift the ZooScan gently and maintain it (e.g., always hold the tray when you lift it to remove the water or it could fall).  
从样品池里回收样品：轻轻地举起 ZooSCAN 并扶住（例：当你把它举起来倒出水时要一直扶着，否则它可能会掉落）。
3. Clean the tray with a squeeze bottle of freshwater or filtered seawater to avoid contamination between the different samples.  
用装淡水或者过滤海水的洗瓶清洗样品池，避免不同样品之间互相污染。

Prepare another sample or clean and DRY the scanning tray using fresh water if you end a scanning session.

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## 9. CONVERT & PROCESS Images and organisms in batch mode

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The image process can be performed just after the scan if you use a powerful computer or during night or lunch time.

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**It is very important to process the scanned images as soon as possible after their acquisition or at least every day. The process permits to rapidly check (see below) that the background and the samples have been correctly acquired.**

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1. Start Zooprocess if necessary, click once on the Z icon and select the tool “CONVERT & PROCESS Images and organisms in batch mode” and then “Convert AND Process Image AND particles (images in RAW folder)”. It is recommended to keep the default settings for the process.  
如果有需要的话可以打开 Zooprocess，点击“Z”键，并且选择“CONVERT & PROCESS Images and organisms in batch mode”工具，然后选择“Convert AND Process Image AND particles (images in RAW folder)”。对于这个过程来说，推荐保持系统的默认设置。
2. At the end of the process, a Zooprocess window appears with a “NORMAL END...” message.  
在处理完之后，一个 Zooprocess 窗口会提示“NORMAL END...”信息。

FOR YOUR INFORMATION : 供参考的信息:

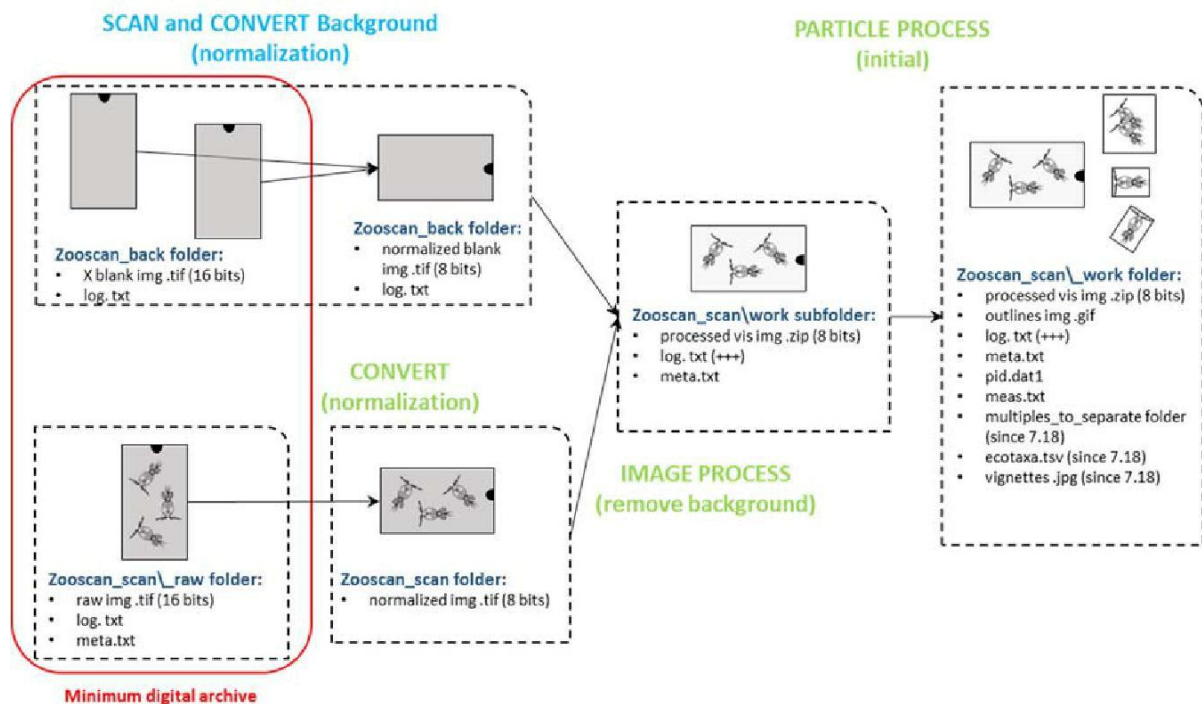
Figure 4 : Description of the standard initial background and image processes. The operations CONVERT, IMAGE PROCESS and PARTICLE PROCESS are all included in the “Convert AND Process Image AND particles (images in RAW folder) » tool.

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- The resulting files are saved in the `_work` and `Pid_results` folders.

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The `.pid` file is a single file that concatenates the `log.txt`, the `meta.txt`, the processing functions applied and the `meas.txt` (table containing all objects (rows) and their measurements (columns)). The measurements that might be utilized to compute the size for the data analysis are Area, Major and Minor (major and minor axes of an ellipse that have the same area of the object measured). Other measurements correspond to variables of shape and texture used for automatic recognition, and of position in the tray. Note that the `PID` file is automatically copied in the `PID_Results` folder of your project.

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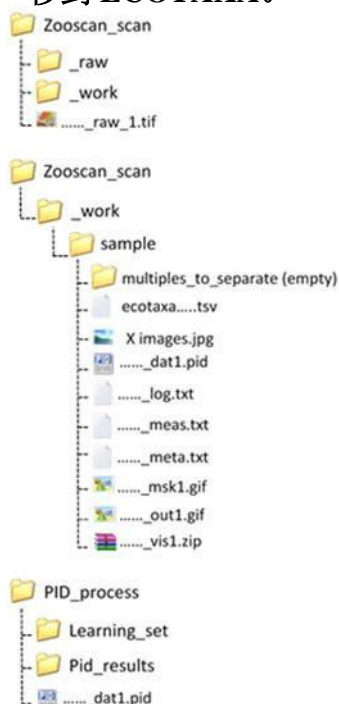
- Since the Zooprocess 7.22 version: vignettes of scanned objects are extracted by default and an “`ecotaxa_*.tsv`” file is created. They are both imported into the ECOTAXA (<http://ecotaxa.obs-vlfr.fr/>) application to predict and classify the organisms. The TSV file contains all data and all useful metadata from the `PID` file (see below).

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- The resulting files remain compatible with the old Plankton Identifier and XnView tools but we **DO NOT** recommend to continue to use them. You can contact us in order to get assistance to move your already classified samples/projects into ECOTAXA.

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## 10. CHECK process by viewing segmented images

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**After any image process**, you must check that the background was well subtracted from the sample image using the dedicated checking tool: “CHECK process by viewing segmented images”. The opened image (“sample\_msk1.gif”) shows if the background was properly extracted from your image, (i.e., no saturated areas, with many dots). You can also check on this image the degree of aggregated organisms.

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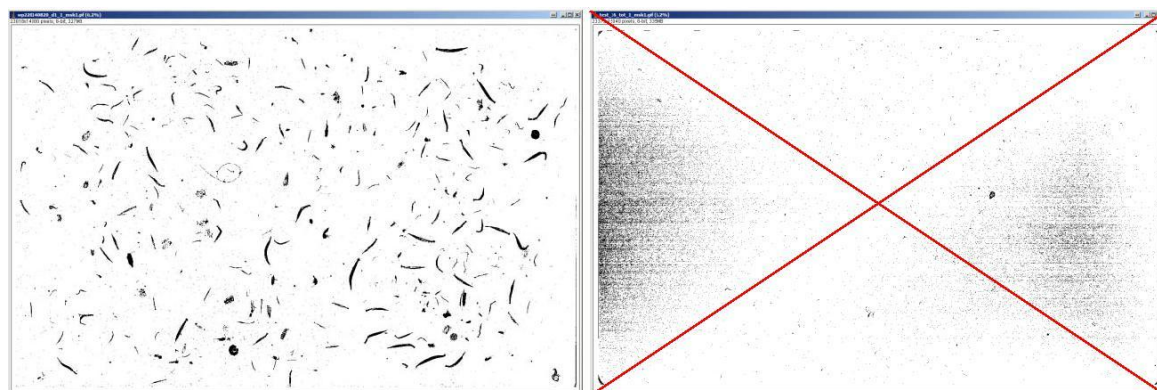
Figure 5 : Specimen of image "masks" : correct and incorrect  
图5：样品图片：正确和错误示范

Some basic reasons for bad quality images can be :

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- Too many organisms  
组织样品过多
- Too many very large organisms or badly positioned  
大型样品太多或者放置位置错误
- Bad background (dirty, condensation, insufficient quantity of water) => all samples of the same day (processed using the same background) present similar defaults  
背景很差（脏、冷凝、水量不够）=> 同一天的所有样品（使用相同的背景进行处理）都存在相同的默认值。
- Bad sample (dirty, condensation, insufficient quantity of water)  
样品质量差（脏、冷凝、水量不够）
- Insufficient delay between the Preview of the sample and its scan  
样品的预览和扫描之间的时间差不足。

## 11. CHECKING “Multiples” and SEPARATION using selected vignettes of multiples 检查“多个对象”并对选定的含有多个对象的小图进行分离

If you are not satisfied with the manual separation that you performed on the scanning tray (too many objects touching in the image) and you notice that too many vignettes contains touching objects, you should perform an additional separation on the vignettes.

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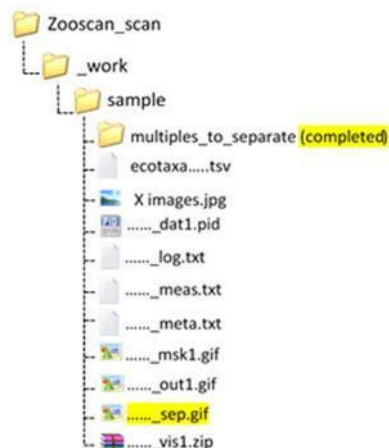
1. Use XnView or any image browser application to check the sample folder (sub folder of the \_work folder) for vignettes containing touching objects and move (cut or copy) them into the “multiples\_to\_separate” subfolder.

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2. Use the tool “SEPARATION using selected vignettes of multiples” to separate the touching organisms by drawing lines between them. The latest version of the tool (Zooprocess 7.25) permits to draw many lines on a single vignette in a batch.

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## 12. RE-PROCESS after separation 分离之后再处理

You must redo the “CONVERT & PROCESS Images and organisms in batch mode” and select “Process again particles from processed images to include SEPARATION MASK (images in WORK sub-folders)” of samples to take benefit of the separation and get results for the resulting objects. At the end of the sample processing a zooprocess window appears with “NORMAL END”.

你必须重新进行“CONVERT & PROCESS Images and organisms in batch mode”步骤，并选择“Process again particles from processed images to include SEPARATION MASK (images in WORK sub-folders)”选项，以利于样品的分离并获得最终的结果。在样品处理的最后，Zooprocess 软件会弹出“NORMAL END”窗口。


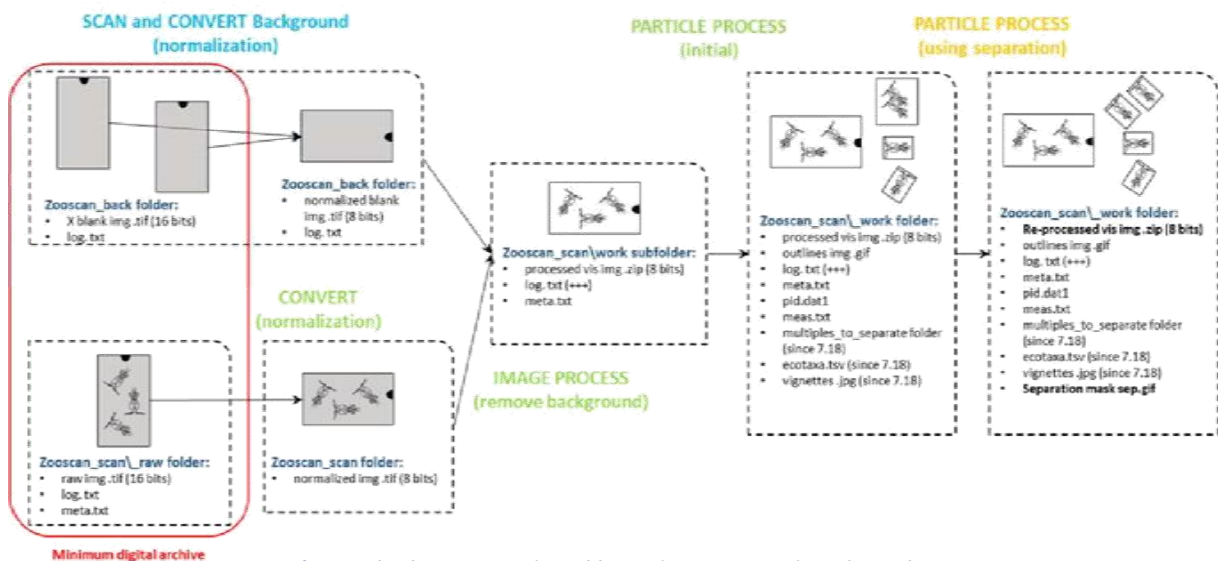
Figure 6 : Standard processes plus additional separation of touching objects  
图6：标准处理程序加互相接触对象额外的分离

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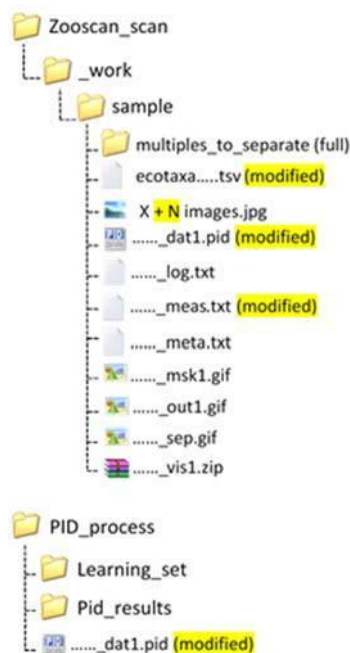The USER mode is set as default when you create a project. It simplifies the daily work limiting the options the user can access, and preventing most of possible manual errors.

After entering a project, you can switch to the ADVANCED mode to get access to the configuration tools and all other options selecting the “SWITCH Mode” tool, wich is at the bottom of the options list.

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